

Amendment for SIRV-Set 3 Lot No. 001602

SIRV108 removed from SIRV isoform mix E0 / ERCC in Lot No. 001602; for compensation SIRV203 is 2x concentrated

Dear SIRV customer,

As part of the SIRV production quality assurance we had to exclude SIRV108 from mixes with the lot numbers 001602 (see labels on the vials). These lots now contain 68 SIRV transcript variants each. For compensation, SIRV203 of similar length and from the same Submix 2 is present in 2-fold concentration to maintain the overall mass and molarity. Its concentration is now 2-fold when compared to Submix 2 and must be treated as a separate entity.

Implications

- Use the annotations as given by [SIRV_Set3_Lot001602_Sequences_170612a \(ZIP\)](#), with the concentration values as shown in [SIRV_Set3_Lot001602_sequence-design-overview-170612a \(XLSX\)](#).
- Exclude SIRV108 from any customized evaluation pipelines, increase the concentration of SIRV203 by a factor of 2, and treat SIRV203 as separate entity (e.g., Submix2.1) for all Submix-related evaluation parameters.


All other expected absolute and relative concentration values are not affected.

File updates

The reference files are accessible from download section, <https://www.lexogen.com/sirvs/downloads/>, by selecting the appropriate SIRV-Set and lot number.

We apologize for any inconvenience this causes. If you have any questions, please contact us at info@lexogen.com or call Lexogen Support at +43(0) 1 345 1212-41.

Sincerely,

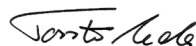

Torsten Reda
Chief Scientific Officer